

# CARSON<sup>®</sup>

MM-450



## POCKETMICRO<sup>™</sup>

### 20x - 60x ZOOM POCKET MICROSCOPE

#### INSTRUCTIONS FOR USE




#### Product Overview

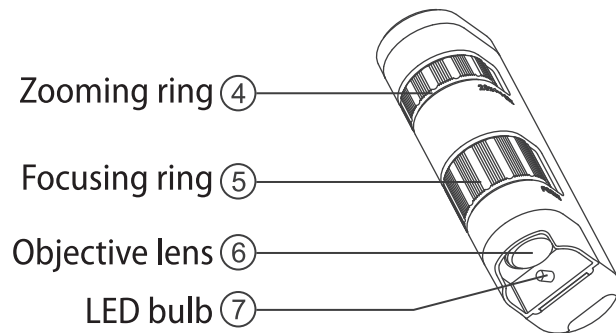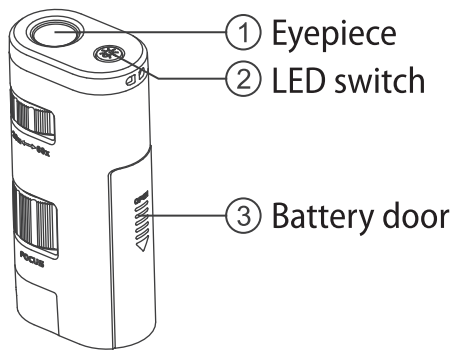


#### Installing the Battery


1. Slide open the battery door ③ as shown below.

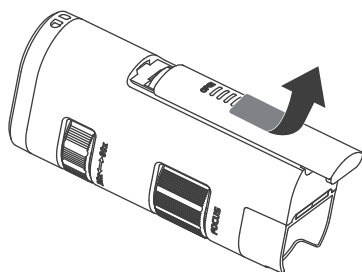

2. Install 1 AA battery with the polarity shown below.

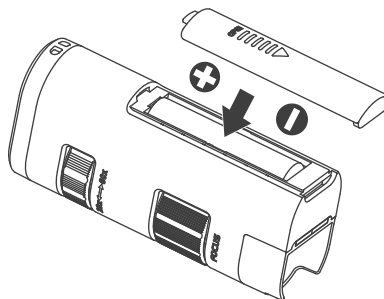


3. Reattach battery door ③ as shown below.

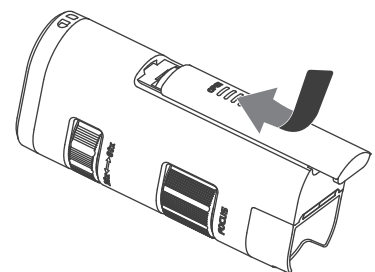


## Using the Microscope


Place the viewing subject directly below the objective lens ⑥.

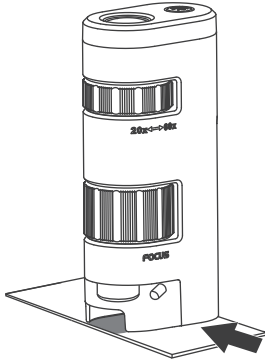


Press the LED button ② to turn on the LED.

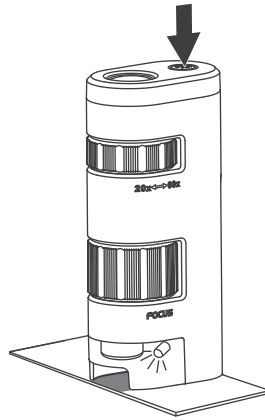


Adjust the zooming ring ④ to the desired level of magnification.

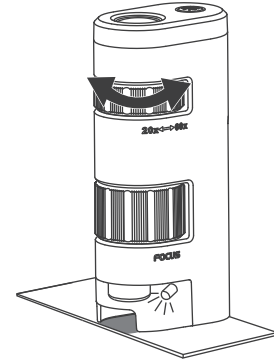

Rotate the focusing ring ⑤ until the image is sharp and clear.

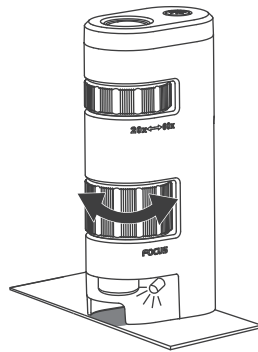

Turn off the LED when finished by pressing the LED button ②.

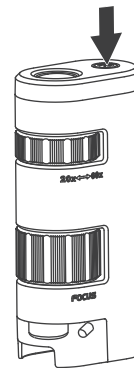

## Specifications

**Magnification:** 20 - 60x  
**Eye Relief:** 8.1mm - 9.7mm  
**Optical Zoom:** 3x  
**Battery Type:** 1 x AA battery (not included)

**Illumination:** LED  
**Dimensions:** 3.9" x 1.8" x 1.1"  
**Weight:** 1.9 oz. (excluding battery)

## Safety Information

1. To avoid possible permanent eye damage, do not look at the sun or any bright objects with the microscope.
2. To avoid possible permanent eye damage, do not place the lighted microscope directly upon human eye.
3. Remove the battery before storing the microscope for an extended period.
4. Inspect all batteries for signs of corrosion or leakage. Do not expose to fire or excessive heat. Use alkaline batteries only, rechargeable batteries not recommended. Insert battery in the direction indicated.
5. Remove exhausted batteries from device.

Öä d ä ~ ö ä / ä ~ K



100 Lauman Lane, Suite A, Hicksville, NY 11801  
Tel: (877) 877-7274 | Fax: (516) 801-2046  
Email: [Info@nyscopes.com](mailto:Info@nyscopes.com)  
[www.microscopeinternational.com](http://www.microscopeinternational.com)