



Meiji Techno America, Inc.

5895 Rue Ferrari
San Jose, CA 95138
Tel: 408.226.3454 | Fax: 408.226.0900
www.meijitechno.com


Öä dā~ cāÄ~K



100 Lauman Lane, Suite A, Hicksville, NY 11801
Tel: (877) 877-7274 | Fax: (516) 801-2046
Email: Info@nyscopes.com
www.microscopeinternational.com

MG800 and MG900 Magnifier Series Instruction Manual

Basic Operations


To turn the unit on, please press the power  button and release.

Special Features

White Light Operation

Press Sequence button to toggle through lighting options:

- Push once, turns on one LED light display
- Second time, turns on two LED light display
- Third time, turns on all three LED light display
- Fourth time, turns off all LED light displays

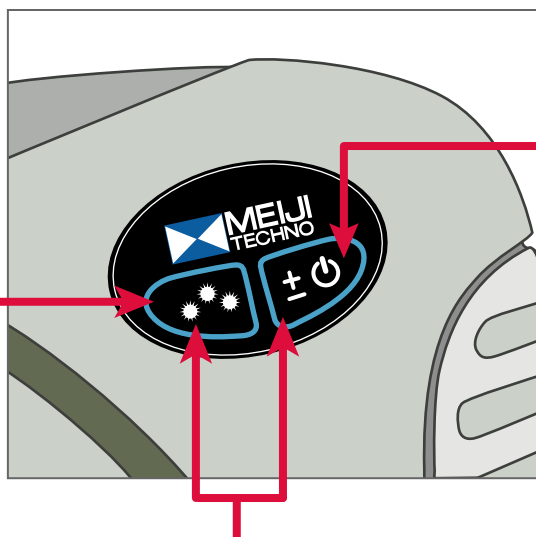

Adjustable Intensity Control

Press and hold Power button down to adjust brightness cycles from high to low and back, continuously. Release button when desired brightness is achieved.

Automatic Timer Function for Light Efficiency

This unit is equipped with an automatic timer function which turns the unit off at 4 hours and 5 minutes, for energy management. To enable timer, press and hold "Both" buttons for 2 seconds and the unit will "Blink once" to acknowledge. To disable the timer, press both buttons, holding again for 2 seconds and the unit will "Blink Twice" to acknowledge.

User setting will be saved until changed

Questions and Concerns?

We at Meiji Techno America carry well designed Green-Lite LED Magnifiers to provide you with years of flawless performance. Should you have questions about these lights, we would love to help you! Further information is available online at meijitechno.com, or you can call us at 408-226-3454, or email us at info@meijitechno.com.

Ask about your different color options:

