OPTIKAM B0.5/B2

# INSTRUCTION MANUAL

Model
4083.B05
4083.B2

Version: 4  
Issued: 28, 04, 2014

---

## Installation of camera drivers and software

Operating system requirements: Windows XP, Windows Vista, Windows 7 or Windows 8. If your PC is Windows XP: before the installation of your Camera please check if Service Pack2 is installed. If not, please update your XP from the Microsoft web site. If your PC has a USB port less than 2.0, there may be some problems in capturing image from the camera due to the slow transmission.

1. Please insert your CD and run:  
    \optika vision lite \ setup.exe
2. Now Optika Vision Lite icon appears on your desktop.
3. Please run \ optika MiPro \ setup.exe
4. Now optika MiPro icon appears on your desktop.
5. Choose from the folder “drivers” the model of camera you want to use.
6. As you can see you don't need to instal any drivers.
7. At the end of the installation process restart the PC if requested.
8. Plug the Camera in your PC: the camera will be identified automatically.

### Some notes

You should know that your Windows PC needs to install a Driver before capturing images from the camera: after installing the driver from the CD, Windows needs to recognize the camera before using it.

Often a PC has some ports on the front and some on the back: the ports on the back are usually faster. We suggest to use these ones.

If you connect the camera into another USB port, Windows needs to recognize its Driver again even if it has already done this operation for the first USB port.

Generally, every time you use a peripheral device and after that you connect your Camera, then Windows needs to recognize the Driver again.

It may happen sometimes that Windows doesn't recognize a Camera using one of your PC USB ports, while gives no problem using another port.

To prevent any installation problem we have put on the CD all the software and drivers in specific folders with an appropriate names.

---

## How to test your Camera

In order to check if your Camera works correctly, it's not necessary to put it into the microscope: it's enough to connect it into the USB port, run Vision Lite using the icon on your desktop, select the right driver and check whether the shown image is not totally black.

### Select the right drivers

It's very easy:

For Optikam B0.5 Camera you have to select the driver:  
USB CAMERA

For Optikam B2 Camera you have to select the driver:  
USB CAMERA

To use Optika Vision Lite you have to tell the Software which is the Driver by selecting the microscope-shaped icon.

To use Optika MiPRO you have to select "capture" and then select the camera from the top right screen list.

### Getting started with Optika Vision Lite, a very easy measuring and documentation software

Run Optika Vision Lite by clicking on its icon on your desktop.

Select the right Driver for capturing the image. Note that after choosing the driver and before capturing the image you can set the brightness, saturation, gamma ecc..  
Then you can capture the image.

The image is now ready for being manipulated using Optika Vision Lite.

### Getting started with optika MiPRO, a complete software for measuring, adjusting, capturing images and video clip.

Run Optika MiPRO by clicking on its icon on your desktop.

Select and choose the right Driver for the camera. Now you can capture a preview of a still image or capture a video.

Note that you can set the brightness, saturation, gamma ecc...and your Live image will change according to the modified values.

The image and video are now ready for being saved.

---

## Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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## OPTIKA® S.r.l.