SZM-SMD

INSTRUCTION MANUAL

MANUEL D'INSTRUCTIONS

Description

The SZM-SMD attachment has been designed specifically for the inspection of solder connections, rework and repair of electronic circuits and for mounting small electronic components. One-touch sliding movement allows the user to quickly change from vertical to oblique viewing.

The attachment is composed of three different parts: A, B, C (see Fig. 1).

Fig. 1

Specifications

Rotatable:	360°
Vertical working distance:	103 mm
Oblique working distance:	103 mm
Oblique angle:	adjustable from 30° to 60°

Assembling of the SZM-ZMD attachment

1. Attach part A (see Fig.1) to the inferior part of the SZM head (see Fig.2).
2. Put the head in the stand of the base.
3. Connect part C to part A of the head.
4. Connect part B to part C using the side screw.
5. The final setup should be similar to the one in Fig. 3.

Fig. 2 The SZM stereo head

Fig. 3 The final setup

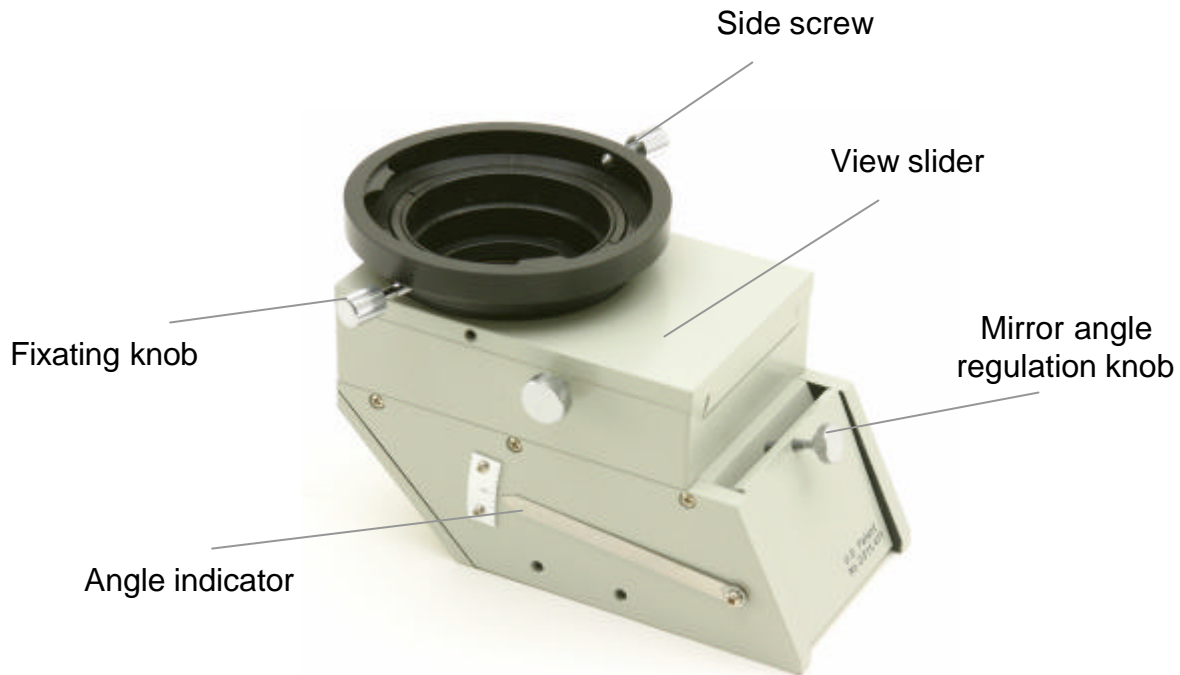
Fig. 4

Switching between oblique and vertical observation

To switch between oblique and vertical view field, loosen the knob on the side of the view slider (see Fig.4). Move the slider horizontally to the view that you prefer and fix it again with the knob.

Regulating the angle of the mirror

To regulate the angle of observation during oblique viewing, turn the mirror angle regulation knob. The angle indicator points at 0 when the mirror is inclined 45°, in its upper position when the inclination is 60° and in the lower position when the mirror is inclined 30°.

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New York Microscope Company Inc